

Verwendung von Histologi@ (digitale Präparate zum Kursus der Mikroskopischen Anatomie an der LMU München) in der Cloud

Univ.Prof. Dr.med. Christoph Schmitz
Inhaber des
Lehrstuhls II für Anatomie
Anatomische Anstalt
Ludwig-Maximilians-Universität München

Laden Sie den kostenlosen Biolucida Viewer unter

<https://biolucida.net>

herunter und installieren Sie ihn auf Ihrem PC oder Mac.

Biolumida Cloud Viewer

Slide Tools Help

Browser

Connected to BiolumidaDemo IC95-FPU-27H1 as BiolumidaDemo

Home

Connections

Tracked Directory:

Showing 1 - 4 of 4 items.

Thumbnail	Name	Description	Tags	Date Modified
	g_1_40.jp2			2013-03-15 14:22:06
	P43_115_GFAP_A06_S0008_2.jp2			2013-03-15 14:22:09
	PS_1_H011_S0014_2.jp2			2013-03-15 14:22:10
	Peruzzi_ImprNett_BH_10302003.jp2			2016-01-26 15:47:06

Showing 1 - 4 of 4 items.

Powered by Biolumida © Copyright 2020 MBP Bioscience, all rights reserved.

Klicken Sie auf „Connections“

Schritt 2: Verbindung mit dem LRZ-Server

Biolumida Cloud Viewer

Slide Tools Help

Browser

Connected to BiolumidaDemo [C95-FPU-27H] as BiolumidaDemo

Home Account

biolumida

Tracked Directories

- ALL
- Classrooms
- 02_Vessels
- ANNOTATED HISTOLOGY ATLAS
- Automatic 3D cell detection
- 3D Virtual Slidebox
- Slide
- Sample Images

Tracked Directory:

Showing 1 - 4 of 4 items

Thumbnail	Name	Description	Tags	Date Modified
	g_1_40.jp2			2013-03-15 14:22:56
	F43_115_GFAP_A06_Slide02.jp2			2013-03-15 14:22:09
	PS_1_H011_Slide14_2.jp2			2013-03-15 14:22:10
	Perazzi_ImprNet_BH_10302003.jp2			2016-01-26 15:47:56

Showing 1 - 4 of 4 items

Biolumida Conn... ? X

QuickCode

Use a QuickCode...

Connections

- BiolumidaDemo
- My Cloud

New Edit Delete

Connect

Featured Sites

- Image Virtual Slidebox
- SPARC Image Collection
- GENEATranscripts Project
- Biolumida Demo Collection

Falls hier noch eine andere Verbindung angezeigt wird (z.B. aus einer früheren Installation), entfernen Sie sie durch Klicken auf „Delete“

Schritt 2: Verbindung mit dem LRZ-Server

Biolucida Cloud Viewer

Slide Tools Help

Browser

Connected to BiolucidaDemo IC95-FPU-27H1 as BiolucidaDemo

Home Account

biolucida

Tracked Directories

- ALL
- Classrooms
- 02_Vessels
- ANNOTATED HISTOLOGY ATLAS
- Automatic 3D cell detection
- 3D Virtual Slidebox
- Image
- Sample Images

Tracked Directory:

Showing 1 - 4 of 4 items.

Thumbnail	Name	Description	Tags	Date Modified
	g_1_40.jp2			2013-03-15 14:22:56
	F43_115_GFAP_A06_Slide02_2.jp2			2013-03-15 14:22:00
	PS_1_H011_Slide14_2.jp2			2013-03-15 14:22:10
	Pericardial_injection_10102003.jp2			2010-01-26 15:47:56

Showing 1 - 4 of 4 items.

Biolucida Conn... ? X

QuickCode

Use a QuickCode...

Connections

BiolucidaDemo

My Cloud

New Edit Delete

Connect

Featured Sites

- Image Virtual Slidebox
- SPARC Image Collection
- GENEACollections Project
- Biolucida Demo Collection

Klicken Sie auf „New“

Biolumida Cloud Viewer

Slide Tools Help

Browser

Connected to BiolumidaDemo [C95-FPU-27H] as BiolumidaDemo

biolumida

Tracked Directories

Tracked Directory:

Thumbnail	Name	Description	Tags	Date Modified
	g_1_40.jpg			2013-03-15 14:22:06
	F43_115_GFAP_A06_SlowR2_2i.jpg			2013-03-15 14:22:09
	PS_1_H011_Sden14_2i.jpg			2013-03-15 14:22:10
	Peruzzi_Ingriffel_BH_10302003.jpg			2016-01-26 15:47:06

Geben Sie die folgenden Informationen ein:

- Connection Name: LRZ
- URL: vm-138-246-234-196.cloud.mwn.de
- Username: Student
- Password: Histokurs2021!

Biolumida Connections

QuickCode

Use a QuickCode...

Connections

biolumidaDemo

My Cloud

Connection Info

Connection Name: LRZ

URL: vm-138-246-234-196.cloud.mwn.de

☐ Connect as guest user

Username: Student

Password: *****

Verify password: *****

Save Cancel

Biolumica Cloud Viewer

Slide Tools Help

Browser

Connected to BiolumicaDemo [C95-FPU-27H] as BiolumicaDemo

biolumica

Tracked Directories

- ALL
- Classrooms
- 02_Vessels
- ANNOTATED HISTOLOGY ATLAS
- Automatic 3D cell detection
- DOWN VIRTUAL SLIDEBOX
- Slide
- Sample Images

Tracked Directory:

Showing 1 - 4 of 4 items.

Thumbnail	Name	Description	Tags	Date Modified
	g_1_40.jp2			2013-03-15 14:22:06
	F4_3_15_GFAP_A06_Slide02_2.jp2			2013-03-15 14:22:09
	PS_1_H011_Slide14_2.jp2			2013-03-15 14:22:10
	Peruzzi_Ingriftell_BH_10302003.jp2			2016-01-26 15:47:06

Showing 1 - 4 of 4 items.

Powered by Biolumica © Copyright 2020 MBP Bioscience, all rights reserved

Biolumica Connections

QuickCode

Use a QuickCode...

Connections

BiolumicaDemo

My Cloud

Connection Info

Connection Name:

LRZ

URL:

vm-138-246-234-196.cloud.mwn.de

☐ Connect as guest user

Username:

Student

Password:

Verify password:

Save Cancel

New Edit Delete

Connect

Abschließend klicken Sie auf „Connect“

Schritt 2: Verbindung mit dem LRZ-Server

Biolumida Cloud Viewer

Slide Tools Help

Browser

Connected to BiolumidaDemo [C95-FPU-27H] as BiolumidaDemo

biolumida

Tracked Directories

- ALL
- Classrooms
- 02_Vessels
- ANNOTATED HISTOLOGY ATLAS
- Automatic 3D cell detection
- DOWN VIRTUAL SLIDEBOX
- Image
- Sample Images

Tracked Directory:

Showing 1 - 4 of 4 items.

Thumbnail	Name	Description	Tags	Date Modified
	g_1_40.jp2			2013-03-15 14:22:06
	F43_115_GFAP_A06_Slide02_2.jp2			2013-03-15 14:22:09
	PS_1_H011_Slide14_2.jp2			2013-03-15 14:22:10
	Peruzzi_Ingriffel_BH_10302003.jp2			2016-01-26 15:47:06

Showing 1 - 4 of 4 items.

Powered by Biolumida © Copyright 2020 MPF Bioscience, all rights reserved

Biolumida Connections

QuickCode

Use a QuickCode...

Connections

BiolumidaDemo

My Cloud

Connection Info

Connection Name:

LRZ

URL:

vm-138-246-234-196.cloud.mwn.de

☐ Connect as guest user

Username:

Student

Password:

Verify password:

Save

Cancel

Connect

Klicken Sie auf „Save“

Biolumida Cloud Viewer

Slide Tools Help

Browser

Connected to BiolumidaDemo [C95-FPU-27H] as BiolumidaDemo

biolumida

Tracked Directories

- ALL
- Classrooms
- 02_Vessels
- ANNOTATED HISTOLOGY ATLAS
- Automatic 3D cell detection
- DOWN VIRTUAL SLIDEBOX
- Image
- Sample Images

Tracked Directory:

Showing 1 - 4 of 4 items.

Thumbnail	Name	Description	Tags	Date Modified
	g_1_40.jp2			2013-03-15 14:22:06
	F4_3_15_GFAP_A06_Slide02_2.jp2			2013-03-15 14:22:09
	PS_1_H011_Slide14_2.jp2			2013-03-15 14:22:10
	Peruzzi_ImprNett_BH_10302003.jp2			2016-01-26 15:47:06

Showing 1 - 4 of 4 items.

Powered by Biolumida © Copyright 2020 MBP Bioscience, all rights reserved

Biolumida Connections

QuickCode

Use a QuickCode...

Connections

- BiolumidaDemo
- LRZ
- My Cloud

Connection Info

Connection Name: LRZ

URL: vm-138-246-234-196.cloud.mwn.de

☐ Connect as guest user

Username: Student

Password: *****

Verify password: *****

Save Cancel

New Edit Delete

Connect

Abschließend klicken Sie auf „LRZ“ und dann auf „Connect“

Hier finden Sie die Präparate des Histokurses, sortiert nach den einzelnen Kurstagen

Biolumida Cloud Viewer

Slide Tools Help

Browser

Connections Connected to LRZ as Student

Home Account

biolumida

Tracked Directories

- all

Classrooms

- 01 Ekt, Die Zelle
- 02 Mikroskopie
- 03 Epithel 1, Epithel
- 04 Epithel 2, Drüsenepithel
- 05 Bindegewebe
- 06 & 07 Knochen, Knochen, Entwicklung, Gelenk
- 08 Muskelgewebe
- 09 Nervengewebe
- 10 Plazenta, Zervix
- 11 Zellorganelle, sekretorischer Weg
- 12 Zahnbildung, Zahnschmelz
- 13 Kreislauforgane
- 14 Blut, HIN, Abwehrzellen
- 15-16 Immunsystem Abwehrorgane
- 17 Auge - visuelles System
- 18 Haut, Ohr
- 19 Atmungsorgane
- 20-22 Endokrine Organe
- 22 Mundhöhle**
- 23 Magen-Darm-Trakt
- 24 Leber, Gallenwege, Pankreas
- 25-26 Niere, Harnwege
- 27-29 Weibliche Geschlechtsorgane, Plazenta, Mamme
- 30 Zahn
- 31 Zahnmedizin
- Sample Images
- 20 Humanmedizin

Classroom: 22 Mundhöhle

Showing 1 - 14 of 14 Items


Select All Medium Thumbnails



Showing 14 - 14 of 14 Items

Powered by Biolumida © Copyright 2020 MHP BioScience, all rights reserved.

Bei Klick auf einen der Kurstage („Classrooms“) zeigt das System die zugehörigen Präparate als Vorschaubilder („thumbnails“). Klicken Sie auf eines dieser Vorschaubilder, wird das entsprechende Präpatat geöffnet.

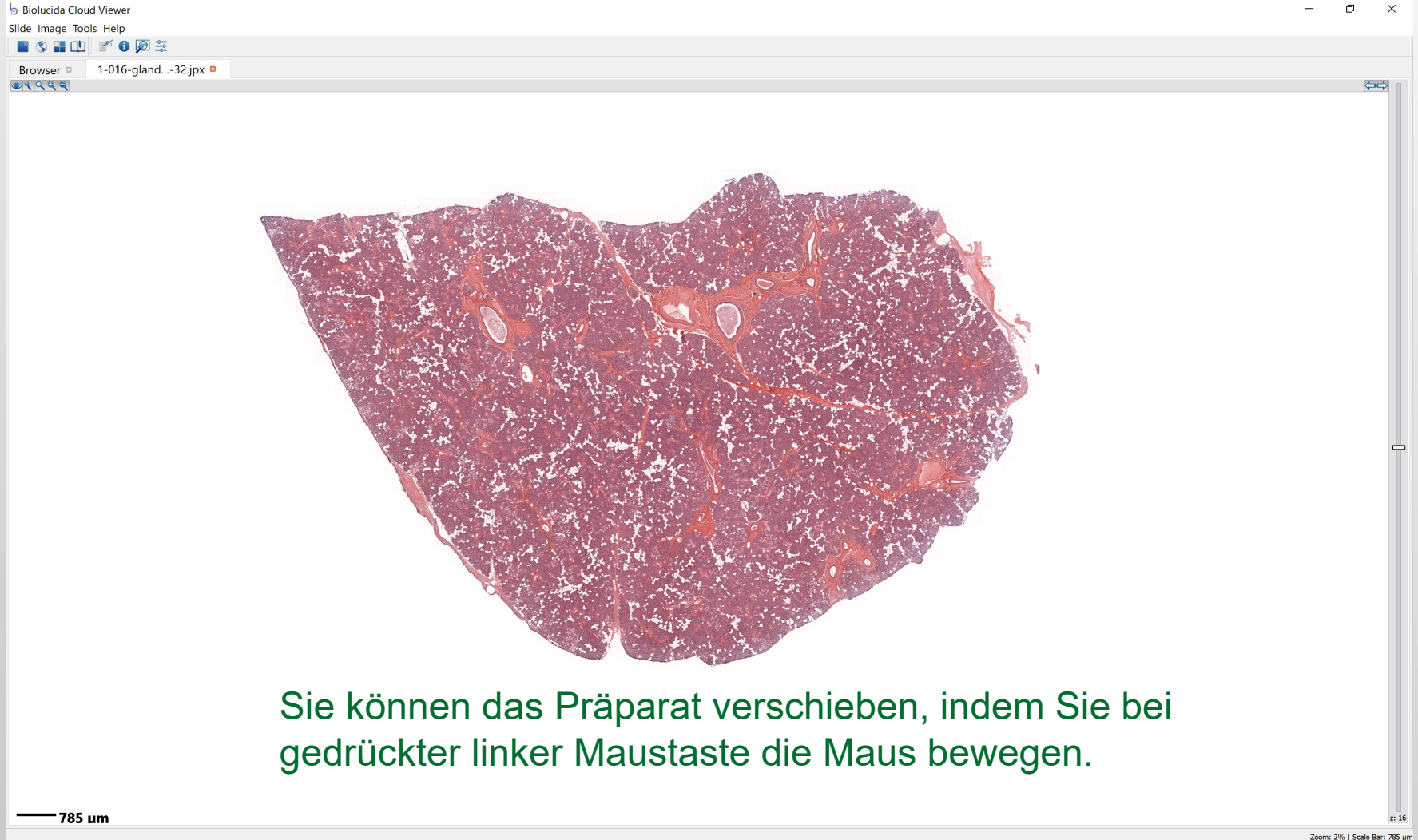



Biolucida Cloud Viewer

Slide Image Tools Help



Browser 1-016-gland...-32.jpg



Zur Beendigung des Studiums eines der Präparate klicken Sie auf das weiße Kreuz in dem roten Kästchen („Close Tab“).

50 µm

Zoom: 25% | Scale Bar: 50 µm